

CE 287

Plate and Frame Filter Press



- * **Separation of solids from suspensions with a plate and frame filter press**
- * **Discontinuous cake filtration**
- * **Practical experiments on a laboratory scale**

concentration of the filtrate to be determined. A drying chamber is recommended for evaluation of the experiments.

The well-structured instructional material sets out the fundamentals and provides a step-by-step guide through the experiments.

Technical Description

Plate and frame filter presses are used in the beverage industry, for example, to clarify intermediate products.

A suspension of diatomite and water (recommended) is prepared in a tank. A pump ensures that the solid remains suspended and does not settle. The pump delivers the suspension into the individual separating chambers of the plate and frame filter press. A separating chamber is formed by one filter frame and two filter plates. The filter plates are grooved and covered over with filter cloths. The filtrate passes through the filter cloth and flows via the grooves in the plates into a collecting pipe. The filtrate exits the plate and frame filter press through the collecting pipe and is collected in the filtrate tank. The solid material is separated off at the filter cloth, where it forms a growing filter cake. As the filter cake becomes thicker, its flow resistance also increases. When the separating chamber is full, or a maximum pressure difference has been reached, the filtration process is ended. The plates and frames of the plate and frame filter press are pulled apart. The filter cake can be removed. For the next filtration the plates and frames must be pushed back together. A spindle is used to press them together. The press forces ensure that the suspension does not leak from the contact points between the plates and the frames, but is forced through the filter cloth.

The flow rate through the plate and frame filter press is adjusted by a valve. The pressure occurring during filtration is indicated on a manometer. The filtrate tank is scaled. This means a stopwatch can be used to measure the flow rate. An included opacimeter allows the solid

Learning Objectives / Experiments


- learning the fundamental principle and method of operation of a plate and frame filter press
- production of a suspension
- removal of the filter cake
- insertion of the filter cloth
- fundamentals of cake filtration: Darcy's equation
- variation in time of filtrate quantity and solid concentration in filtrate
- mass of filter cake dependent on filtrate quantity

CE 287


Plate and Frame Filter Press



1 switch cabinet with controls, 2 suspension tank, 3 filtrate tank outlet and overflow, 4 filtrate tank, 5 spindle, 6 plate and frame filter press

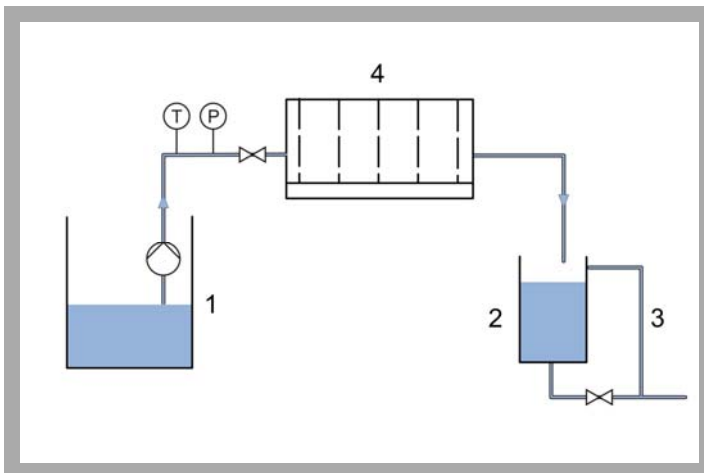

1 tank with pump, 2 filtrate tank, 3 overflow, 4 plate and frame filter press; T thermometer, P manometer

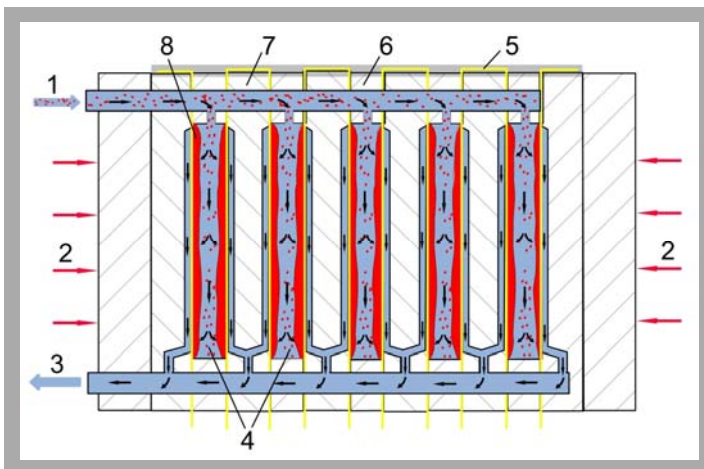

Fundamental principle of a plate and frame filter press: 1 suspension inlet, 2 press forces, 3 filtrate outlet, 4 separating chambers, 5 filter cloth, 6 filter frame, 7 filter plate, 8 filter cake

Specification

- [1] plate and frame filter press for discontinuous cake filtration
- [2] HDPE tank to produce a suspension
- [3] centrifugal pump to deliver the suspension to the plate and frame filter press
- [4] plate and frame filter press with 10 opening separating chambers for removal of the filter cake
- [5] PMMA tank with level scale for filtrate
- [6] adjustment of suspension flow rate by valve
- [7] thermometer and manometer in inlet
- [8] portable opacimeter to measure the solid concentration in the filtrate

Technical Data

- Plate and frame filter press
 - filter area: approx. 0,72m²
 - working pressure: approx. 0,4...2,5bar
- Centrifugal pump (submersible pump)
 - max. flow rate: 4,5m³/h
 - max. head: 45m
- Tanks
 - suspension tank: 200L
 - filtrate: 20L
- Measuring ranges
 - pressure: 0...4bar
 - temperature: 0...60°C

Dimensions and Weight

- LxWxH: 1.900x790x1.900mm
- Weight: approx. 190kg

Connections

- 230V, 50/60Hz, 1 phase or 120V, 60Hz/CSA, 1 phase
- Water connection

Scope of Delivery

- 1 plate and frame filter press
- 1 portable opacimeter
- 1 stopwatch
- 1 set of filter cloths
- 2 hoses
- 1 set of instructional material

Order Details

083.28700 CE 287 Plate and Frame Filter Press